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1. Zhao J, Xu S, Song F, Nian L, Zhou X, Wang S. 2,3,5,4'-tetrahydroxystilbene-2-O-β-D-glucoside protects human umbilical vein endothelial cells against lysophosphatidylcholine-induced apoptosis by upregulating superoxide dismutase and glutathione peroxidase. IUBMB Life. 2014 Oct;66(10):711-22.
2. Zhao H, Wang R, Tao Z, Yan F, Gao L, Liu X, Wang N, Min L, Jia Y, Zhao Y, Ji X, Luo Y. Activation of T-LAK-cell-originated protein kinase-mediated antioxidation protects against focal cerebral ischemia-reperfusion injury. FEBS J. 2014 Oct;281(19):4411-20.
3. Li X, Xu L, Zhou W, Zhao Q, Wang Y. Chronic exposure to microcystin-LR affected mitochondrial DNA maintenance and caused pathological changes of lung tissue in mice. Environ Pollut. 2015 Dec 16;210:48-56.
4. Liu J, Tang X, Wang Y, Zang Y, Zhou B. A Zostera marina manganese superoxide dismutase gene involved in the responses to temperature stress. Gene. 2016 Jan 10;575(2 Pt 3):718-24.
5. Chen W, Su H, Xu Y, Bao T, Zheng X. Protective effect of wild raspberry (Rubus hirsutus Thunb.) extract against acrylamide-induced oxidative damage is potentiated after simulated gastrointestinal digestion. 2016 Apr 1;196:943-52.
6. Lin J, Zou Y, Cao K, Ma C, Chen Z. The impact of heterologous catalase expression and superoxide dismutase overexpression on enhancing the oxidative resistance in Lactobacillus casei. J Ind Microbiol Biotechnol. 2016 May;43(5):703-11.
7. Guo H, Zhang N, Liu D, Wang P, Ma X. Inhibitory effect on the proliferation of human hepatoma induced by cell-permeable manganese

superoxidodismutase. *Biomed Pharmacother.* 2016 Oct;83:1379-1386.

8. Zhuang C, Xu NW, Gao GM, Ni S, Miao KS, Li CK, Wang LM, Xie HG. Polysaccharide from *Angelica sinensis* protects chondrocytes from H₂O₂-induced apoptosis through its antioxidant effects in vitro. *Int J Biol Macromol.* 2016 Jun;87:322-8.
9. Sun X, Jiao X, Ma Y, Liu Y, Zhang L, He Y, Chen Y. Trimethylamine N-oxide induces inflammation and endothelial dysfunction in human umbilical vein endothelial cells via activating ROS-TXNIP-NLRP3 inflammasome. *Biochem Biophys Res Commun.* 2016 Dec 2;481(1-2):63-70.
10. Sun Y, Xiu C, Liu W, Tao Y, Wang J, Qu YI. Grape seed proanthocyanidin extract protects the retina against early diabetic injury by activating the Nrf2 pathway. *Exp Ther Med.* 2016 Apr;11(4):1253-1258.
11. Pan K, Li X, Chen Y, Zhu D, Li Y, Tao G, Zuo Z. Deferoxamine pre-treatment protects against postoperative cognitive dysfunction of aged rats by depressing microglial activation via ameliorating iron accumulation in hippocampus. *Neuropharmacology.* 2016 Dec;111:180-194.
12. Yang J, Liu X, Zhang X, Jin Q, Li J. Phenolic Profiles, Antioxidant Activities, and Neuroprotective Properties of Mulberry (*Morus atropurpurea* Roxb.) Fruit Extracts from Different Ripening Stages. *J Food Sci.* 2016 Oct;81(10):C2439-C2446.
13. Wang J, Chen X, Wang F, Zhang J, Li P, Li Z, Xu J, Gao F, Jin C, Tian H, Zhang J, Li W, Lu L, Xu GT. OFD1, as a Ciliary Protein, Exhibits Neuroprotective Function in Photoreceptor Degeneration Models. *PLoS One.* 2016 May 19;11(5):e0155860.
14. Cao XN, Shen LJ, Wu SD, Yan C, Zhou Y, Xiong G, Wang YC, Liu Y, Liu B, Tang XL, Guo M, Liu DY, Long CL, Sun M, He DW, Lin T, Wei GH. Urban fine particulate matter exposure causes male reproductive injury through destroying blood-testis barrier (BTB) integrity. *Toxicol Lett.* 2017 Jan 15;266:1-12.
15. Huang W, Guo HL, Deng X, Zhu TT, Xiong JF, Xu YH, Xu Y. Short-Chain Fatty Acids Inhibit Oxidative Stress and Inflammation in Mesangial Cells Induced by High Glucose and Lipopolysaccharide. *Exp Clin Endocrinol Diabetes.* 2017 Jan 3. doi: 10.1055/s-0042-121493. [Epub ahead of print]
16. Chen W, Su H, Xu Y, Jin C. In vitro gastrointestinal digestion promotes the protective effect of blackberry extract against acrylamide-induced oxidative stress. *Sci Rep.* 2017 Jan 13;7:40514.
17. Hu W, Wang H, Liu Z, Liu Y, Wang R, Luo X, Huang Y. Neuroprotective effects of lycopene in spinal cord injury in rats via antioxidative and anti-apoptotic pathway. *Neurosci Lett.* 2017 Feb 3;642:107-112.

Version 2018.03.08